

FULL SCALE CALIBRATION AND ZERO ERROR RECORD
(Analytical and Precision Weighing Balance)

Unit : X

Department : QC

Balance Code No. : BAL-006

Make : METTLER TOLEDO

Std. Weight Box ID. : WB-002

Month and year : October 2025

Capacity (a) (g) : 300.000

Model : MS303S

Least Count (e) (g) 0.001

Std. Weight Box Certificate No. 24-07-2025-002

Std. Weight Box Valid Upto 2039-12-31

Tolerance: +/- Least count of the balance OR +/- 0.1 % of Standard weight whichever is higher.

Standard weights for calibration (g)		Tolerance (g)	Accuracy Limits(Std wt +/- Tolerance) (g)
Lower(20.e)/(5% of a)/Minimum capacity of the checkweigher / weighing balance	10.000	± 2.000	8.000 - 12.000
Middle-I(20% of a)	20.000	- 3.000 + 2.000	17.000 - 22.000
Middle-II(50% of a)	30.000	- 2.000 + 3.000	28.000 - 33.000
Upper(80% of a)	40.000	- 2.000 + 3.000	38.000 - 43.000

- Note:**
1. Adjust zero error before starting the weighing operation.
 2. Check against standard weights daily before starting the activity.
 3. If more than one standard weight is used for verification then record the standard weight / identification No. of all individual weights used for verification. (E.g. For verification of 200g, if three weights viz. 50g, 50g & 100g is used then record standard weight/identification No. of all the three weights)

Date	Time	Spirit Level (Ok/ Not Ok)	Zero Error	Standard weights kept on the balance (g)		Readings shown (g)	Remark
10.10.2025	11:11	OK	Nil	Lower(20.e)/(5% of a)/Minimum capacity of the checkweigher / weighing balance	10.000	11.203	Ok
				Middle-I(20% of a)	20.000	20.260	Ok
				Middle-II(50% of a)	30.000	30.225	Ok
				Upper(80% of a)	40.000	40.203	Ok

FULL SCALE CALIBRATION AND ZERO ERROR RECORD
(Analytical and Precision Weighing Balance)

Cipla
Goa , Unit X

Preview
Only

Preview
Only

FULL SCALE CALIBRATION AND ZERO ERROR RECORD
(Analytical and Precision Weighing Balance)

2. Uncertainty test:

Sr. No.	Standard Weight (g)	Reading displayed on the balance (g)	Remarks
1	10.000	11.235	Ok
2	10.000	11.235	Ok
3	10.000	11.235	Ok
4	10.000	11.235	Ok
5	10.000	11.235	Ok
6	10.000	11.235	Ok
7	10.000	11.235	Ok
8	10.000	11.235	Ok
9	10.000	11.235	Ok
10	10.000	11.235	Ok
Mean		11.235	-
Standard Deviation (S.D.)		0.0000	-

Uncertainty =
$$\frac{3 \times \text{S.D.}}{\text{Standard Weight (g)}} = 0.000$$

Acceptance criteria : Value of uncertainty should be NMT 0.001 g.

Remark : Complies

FULL SCALE CALIBRATION AND ZERO ERROR RECORD
(Analytical and Precision Weighing Balance)

3. Repeatability test:

Sr. No.	Standard Weight (g)	Reading displayed on the balance (g)	Remarks
1	10.000	11.239	Ok
2	10.000	11.239	Ok
3	10.000	11.239	Ok
4	10.000	11.239	Ok
5	10.000	11.239	Ok
6	10.000	11.239	Ok
7	10.000	11.239	Ok
8	10.000	11.239	Ok
9	10.000	11.239	Ok
10	10.000	11.239	Ok
Mean		11.239	-
Standard Deviation (S.D.)		0.0000	-

Repeatability (%) = $\frac{2 \times (0.41 \times d)}{\text{Standard weight (g)}} \times 100 = 0.01$



Acceptance criteria : Repeatability is NMT 0.10 %

Remark : Complies

FULL SCALE CALIBRATION AND ZERO ERROR RECORD (Analytical and Precision Weighing Balance)

4. Eccentricity test :

Note: Adjust zero error before commencing the weighing operation .

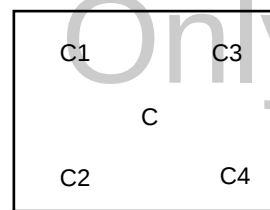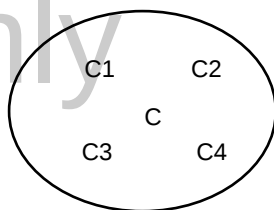

Sr No.	Standard Weight (g)	Position	Reading displayed by the balance (g)	Deviation from center =X-C	Remark
1	10.000	C	11.629	-----	Ok
2	10.000	C1	11.629	0.000	Ok
3	10.000	C2	11.629	0.000	Ok
4	10.000	C3	11.629	0.000	Ok
5	10.000	C4	11.629	0.000	Ok

Where, X=C1, C2, C3, C4 i.e. different corners of the balance.

Acceptance criteria : Deviation is not more than 0.05%.

Remark: Complies

Comment: _____

	Signature	Date & Time
Performed By	Mangesh (11)	10.10.2025 & 11:11:45
Checked By	kuldeep (Z)	10.10.2025 & 11:18:39
Checked By (Quality Assurance)		